

Interpretation guide

The 3M[™] Petrifilm[™] Rapid Yeast and Mold Count Plate is a sample-ready-culture medium system which contains nutrients supplemented with antibiotics, a cold-water-soluble gelling agent and an indicator system that facilitates yeast and mould enumeration.







Yeast vs. mould colonies

To differentiate yeast and mould colonies on the 3M[™] Petrifilm[™] Rapid Yeast and Mold Count Plates, look for one or more of the following characteristics:



Yeast count = 44

The colonies are examples of characteristic **yeast**: small colonies, colonies have defined edges, pink-tan to blue-green in color, colonies appear raised (3 dimensional) and colonies have a uniform colour.



Mould count = 12

The colonies are examples of characteristic **mould**: large colonies, colonies have diffuse edges, blue-green to variable upon prolonged incubation, colonies appear flat and colonies have a dark center with diffused edge.

Growth and colony formation

Incubate Petrifilm Rapid Yeast and Mold count plates at $25^{\circ}C \pm 1^{\circ}C$ or $28^{\circ}C \pm 1^{\circ}C$ for a minimum of $48-60h^{*}$ in a horizontal position with the clear side up in stacks of no more than 40. Certain food types may exhibit clearer growth and colony formation at $28^{\circ}C$.



*If colonies appear faint, allow an additional 12 hours of incubation time for enhanced interpretation. See product instructions for third party validated methods. **The presence of small air bubbles will not prevent accurate counts.

Rapid Yeast and Mold Count Plate

Enzymatic reaction

Food samples may occasionally show interference on the 3M[™] Petrifilm[™] Rapid Yeast and Mold Count Plates, for example:



Count = 0

A uniform blue background colour (often seen from the organisms used in cultured products) should not be counted as TNTC.



Count = 5

Some foods containing high levels of enzymes may cause a uniform blue background. Colony growth will still be visible if an enzyme reaction occurs.



Count = 0 A plate without an enzymatic reaction.



Count = TNTC

Plates containing greater than 150 colonies can either be estimated or recorded as too numerous to count (TNTC).

For a more accurate count, further dilution of sample may be necessary.

Reminders for use

Storage





Store the unopened 3M[™] Petrifilm[™] Rapid Yeast and Mold Count Plate pouches at frozen or refrigerated temperature equal to -20 to 8°C (-4 to 46°F). Use before expiration date on package. It is best to allow pouches to reach room temperature before opening.





Inoculation



Place the Petrifilm Rapid Yeast and Mold count 3 plate on a flat, level surface. Lift the top film and with the pipette perpendicular dispense 1mL of sample suspension onto the center of bottom film.

Roll the top film down onto the sample.

Incubation



Incubate Petrifilm Rapid Yeast and Mold plate at 25°C ± 1°C or 28°C ± 1°C for 48h ± 2h in a horizontal position with the clear side up in stacks of no more than 40. Please refer to the product instructions for third party validated methods.



2

Spreader (6425) or other flat spreader on the center of the Petrifilm Rapid Yeast and Mold count plate.

prototio-Inte



Certain slower growing yeasts and moulds may appear faint. To enhance interpretation of these moulds allow for an additional 12 hours of incubation time.

3M Food Safety offers a full line of products to accomplish a variety of your microbial testing needs. For more product information, visit us at 3M.com/foodsafety/Petrifilm

9



8

3M UK PLC Charnwood Campus 10 Bakewell Road Loughborough LE11 5RB

emeafoodsafety@mmm.com www.3M.co.uk/foodsafety

 		_	
•••	:•. •.•		

6

Press firmly on the center of the

spreader to distribute the sample

evenly. Spread the inoculum over

and Mold count plate growth area

before the gel is formed. Do not slide the spreader across the film.

the entire Petrifilm Rapid Yeast



Use appropriate sterile diluents

to form.

Butterfield's phosphate buffer, buffered peptone water (ISO), 0.1% peptone water, peptone salt diluent, saline solution (0.85-0.90%), bisulphite-free letheen broth or distilled water.

Remove the spreader and leave

count plate undisturbed for at

the Petrifilm Rapid Yeast and Mold

least one minute to permit the gel

Do not use diluents containing citrate, bisulphite or thiosulfate with Petrifilm Rapid Yeast and Mold count plates; they can inhibit growth.

If citrate buffer is indicated in the standard procedure, substitute with 0.1% peptone water, warmed to 40-45°C

User's responsibilities: 3M[™] Petrifilm[™] Plate performance has not been evaluated with all combinations of microbial flora, incubation conditions and food matrices. It is the user's responsibility to determine that any test methods and results meet the user's requirements. Should re-printing of this Interpretation guide be necessary, user's print settings may impact picture and colour quality.

For detailed CAUTIONS, DISCLAIMER OF WARRANTIES/LIMITED REMEDY and LIMITATION OF 3M LIABILITY, STORAGE AND DISPOSAL information and INSTRUCTIONS FOR USE, see Product's package insert.

3M and Petrifilm are trademarks of the 3M company. © 3M 2019. All rights reserved. 70-2011-5050-8. J451005.

Rapid Yeast and Mold Count Plate